


## 2. 제품 사양

MODEL		Magnetic Stirrer			
		MS120D	MS150D	MS200D	MS330D
Stirring	Speed	100 ~ 1500rpm			100 ~ 1300rpm
	Speed display	Digital			
	Capacity(H <sub>2</sub> O)	up to 1L	up to 2L	up to 5L	up to 15L
Speed controller		Feedback control			
Top plate		Ceramic coated top plate			
Dimensions	Plate area	120 x 120	150 x 150	180 x 180	300 x 300
	Overall(WxDxH)	160 x 280 x 107	160 x 280 x 107	200 x 310 x 107	310 x 430 x 120
Electrical supply		AC220V 50/60Hz or AC110V 50/60Hz (option)			
Weight (kg)		2.4	2.5	3.4	5.8

## 3. 각 부의 명칭 및 설명

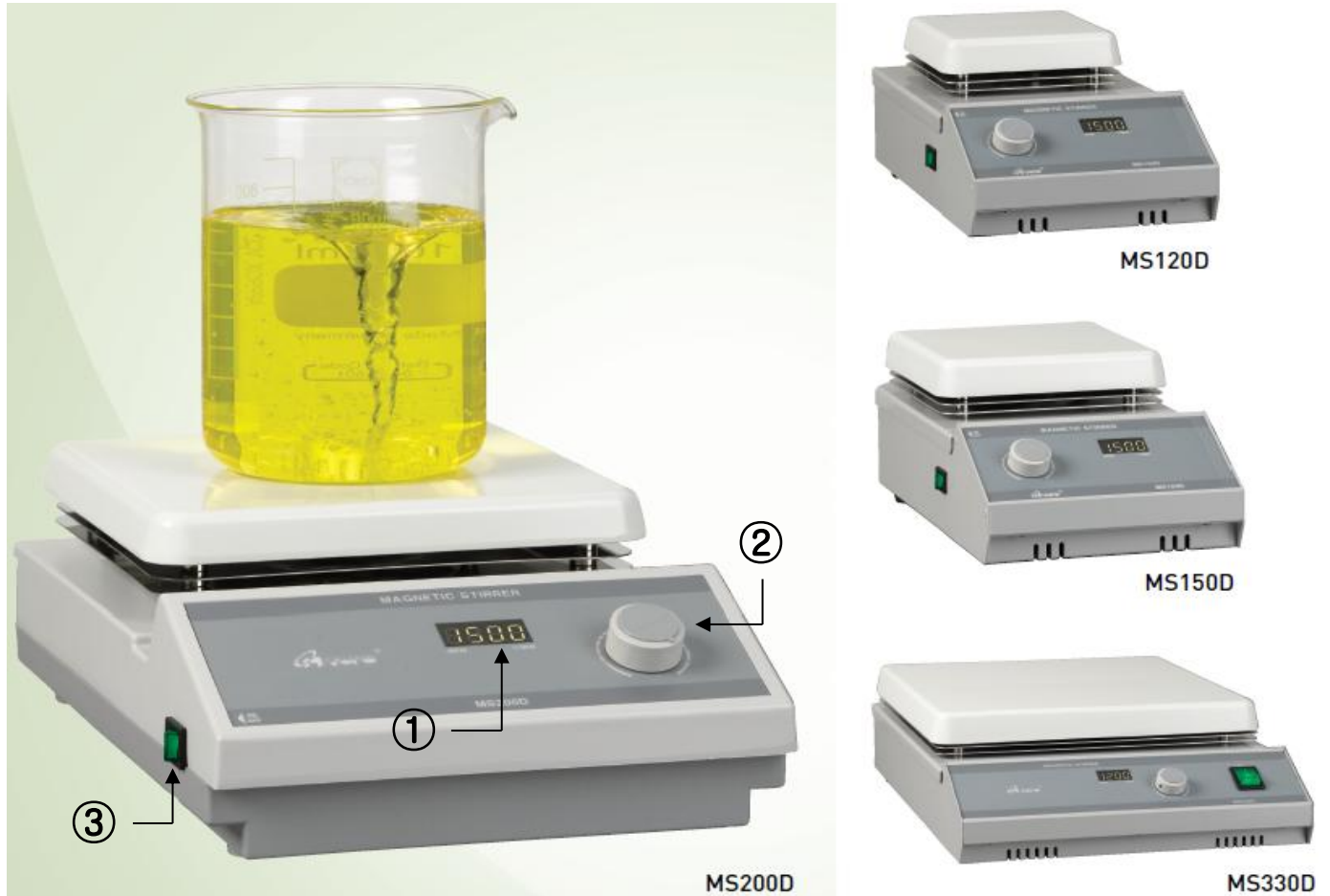

1. Motor 속도 : Motor의 현재rpm과 설정rpm(점멸상태)을 표시합니다.
2. Motor control : Motor 속도를 조절할때 사용합니다.
3. Main power S/W : 주전원을 켜거나 끌때 사용합니다.

## 4. 사용방법

1. 교반:속도볼륨(2)을 돌려서 원하는 속도를 설정하여 사용한다.
2. 실험종료후 실험대상을 제거하고 Main power S/W를 "OFF" 시킨다.
3. 장시간 사용하지 않으실경우 Main power cord를 빼서 청결하게 보관하십시오.