


2. Specifications

MODEL	Electronic overhead stirrer	
	HM1200D	HD1200D
Working capacity	500 ~ 2000ml	500 ~ 5000ml
Motor (out put)	142W	
Speed range	1000 ~ 11000rpm	1000 ~ 12000rpm
Speed controller	Feed-back control	
Speed display	Digital	
Dimensions (WxDxH)	70 x 180 x 170mm	
Weight	3.2kg	2.8kg
Electrical supply	AC220V 50/60Hz	

3. Functional Description

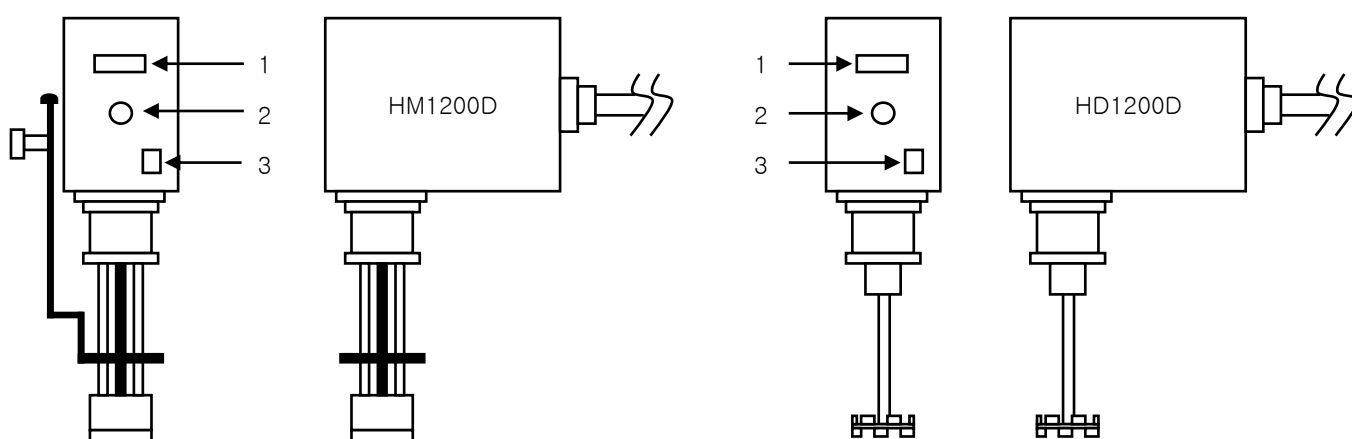

1. Digital display : Display of working RPM and setting value

Turn No.2 dial to see the setting value, then display will show it with blinking

(Actual rpm is 10 times of the displayed value, for example, if the displayed value is 150, actual value is 1500 rpm.)

2.Speed control : Turn the volume to set the speed.

3.Power switch : Use for power ON/OFF. Lamp illuminates on power ON.

4. How to use

1. Assemble the unit.
2. Connect the power cord in rear side to main power supply.
3. Power on with No. 3 switch.
4. Set the required speed turning No.2 volume with looking at No. 2 volume.
5. After finishing a experimentation, place No.2 volume on Min. and separate power cord from the unit.