



2. 각 부의 명칭

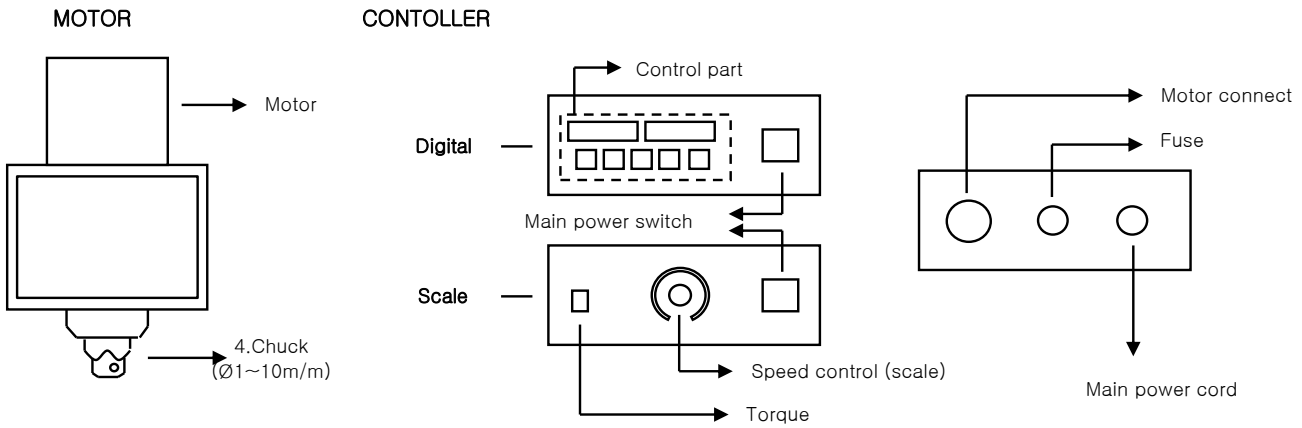

3. 사양

MODEL	Electronic overhead stirrer (BL2000 series)							
	BL2015	BL2015D	BL2010	BL2010D	BL2006	BL2006D	BL2003	BL2003D
Stirring capacity (H2O)	80ℓ		100ℓ		120ℓ		150ℓ	
Viscosity (Max.)	50,000mpas (medium)		100,000mpas (med,high)		150,000mpas (high)		high viscosity	
Motor (out put)	Brushless DC motor 200W							
Torque (Max.)	13kg . cm (127Ncm)		19.5kg . cm (191Ncm)		32kg . cm (318Ncm)		65kg . cm (637Ncm)	
Speed range	65 ~ 1,500rpm		50 ~ 1,000rpm		35 ~ 600rpm		25 ~ 300rpm	
Speed controller	Feed-back control							
Speed display	Scale	Digital	Scale	Digital	Scale	Digital	Scale	Digital
Stirring motion	—	1.Right	—	1.Right	—	1.Right	—	1.Right
		2.left		2.left		2.left		2.left
		3.Auto reverse		3.Auto reverse		3.Auto reverse		3.Auto reverse
Timer	—	99hr 59min	—	99hr 59min	—	99hr 59min	—	99hr 59min
Chuck range	Ø1 ~ 10mm							
Dimensions (WxDxH)	x x							
Weight	kg							
Electrical supply	AC220V 50/60Hz							

4. Scale Type 조작방법

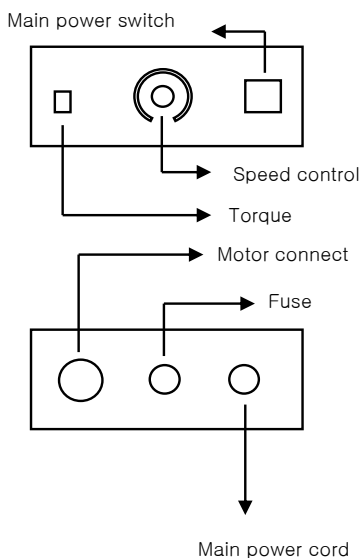

1. 장치를 완성시킨다.
2. 척에 임펠로드를 연결한다.동작도중 떨어지지 않게 단단히 고정한다.
3. 뒤쪽 Motor콘넥터에 Motor본체의 콘넥터를 연결한다.
4. 코오드를 콘센트에 연결한다.
5. 전원스위치를 ON한다.
6. 볼륨을 돌려 원하는 속도를 설정한다.
7. 실험이 끝난후엔 반드시 볼륨을 Min.에 위치시키고 전원코오드를 분리시킨다.

Torque의 표시등이 3개가 깜빡이면 Motor측에 부하가 많이 걸림을 표시합니다.
깜빡거리후 약10초후에 자동정지됩니다.
원인을 찾아 제거후 다시 사용하시기 바랍니다.